

MCS Series PES 0.2µm – Single Layer Filter Capsules

The MCS Series Pharmaceutical Grade Cartridges are designed for product sterility. The allpolypropylene construction provides excellent chemical compatibility and superior flow per unit area as compared to other membrane cartridges. MCS Series Pharmaceutical Grade Cartridges have been designed for final filtration in biopharmaceutical applications. The hydrophilic PES membrane does not require pre-wetting agents, thereby eliminating a potential source of contamination. Also, the PES membrane has low protein binding characteristics to maximize yields. No adhesives, binders, or surfactants are used in the manufacturing process. All cartridges are rinsed with pyrogen-free water to reduce extractables and downtime. All filter cartridges are 100% integrity tested to ensure filter performance and quality. MCS Series Pharmaceutical Grade Cartridges are well-suited for critical applications where superior flow and bacterial retention are required. The filters are manufactured in accordance with GMP and comply with <USP 797> guidelines. The filters are flushed with pharmaceutical-grade purified pyrogens-free water, and 100% integrity tested prior to dispatch. Our filters and free from animal-derived components raw materials can be traced by Lot Number, and are strictly regulated, maintaining ISO 13485 & ISO 14001 quality standards.

MCS Series Filter Cartridge Filters with Code 7: 226 O-Ring with Spear





International Filter Products Revolutionary systems for your industry



International Filter Products

Revolutionary systems for your industry

Application

- Buffers and Media
- LVP (Large Volume Parenteral)
- Vaccines
- Scale up processing

- Product Sterilization
- Bio Bags Compatible
- > Biologics
- Ultrapure Water

Technical Data Sheet

Micron Rating:

Final Membrane: 0.2 µm PES [Sterilizing Grade Membrane]

Effective Filtration Area:

- **5"**; 3.77 *ft*², (0.35 *m*²)
- **10"**; 7.53 ft^2 , (0.7 m^2)
- **20"**; 15.07 *ft*², (1.4 *m*²)
- **30"**; 30.14 *ft*², (2.8 *m*²)
- **40"** ; 60.28 ft^2 , (5.6 m^2)

Materials of Construction:

Membrane: Pharmaceutical Grade PES (Polyethersulfone) Sterilizing Grade

Membrane Feature: Hydrophilic, Low Protein Binding, High Flow Rate

Media Supports: Polypropylene

Shell, Cage, Core, End Caps: Polypropylene

O-Rings: Silicone (Standard)

Sealing: Thermally Bonded

Operating Conditions:

Maximum Forward Differential Pressure:

Liquid: 5.5 bar (80psi) at 77°F/25°C

Gas: 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C

Maximum Operating Temperature: 176°F/80°C




International Filter Products Revolutionary systems for your industry

- Fermentation Broths
- Pharmaceuticals
- > Antibiotics
- Serums




Nominal Dimension:

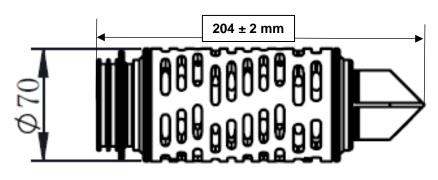

MCS – 5 inches

Part Number	Internal Filter Length (inches)	End Modification	Gasket Material			
MCS2075S	5"	Code 7: 226 O-Ring with Spear	Silicone			
 Find the complete list of available standard inlet and outlet fittings in page 8. ETO- Sterilization is available for all 5-inch cartridges. Add "-ETO" with the original Part Number 						

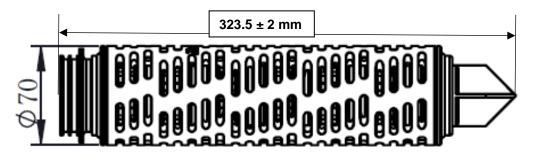

MCS – 10 inches


Part Number	Internal Filter Length (inches)	End Modification	Gasket Material		
MCS2071S	10"	Code 7: 226 O-Ring with Spear	Silicone		
 Find the complete list of available standard inlet and outlet fittings in page 8. ETO- Sterilization is available for all 10-inch cartridges. Add "-ETO" with the original 					

Part Number




International Filter Products

Revolutionary systems for your industry

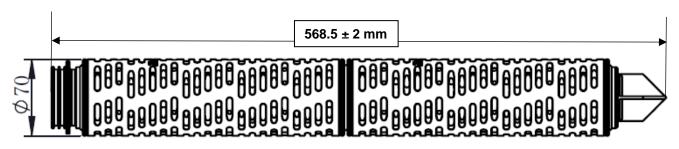

MCS – 20 inches

Part Number	Internal Filter Length	End Modification	Gasket Material	
	(inches)			
MCS2072S 20"		Code 7: 226 O-Ring with Spear	Silicone	
Find the complete list of available standard inlet and outlet fittings in page 8.				

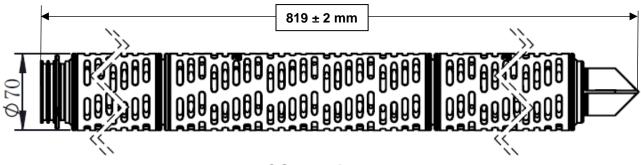


MCS – 30 inches

Part Number	Internal Filter Length	End Modification	Gasket Material		
	(inches)				
MCS2073S	30"	Code 7: 226 O-Ring with Spear	Silicone		
Find the complete list of available standard inlet and outlet fittings in page 8.					





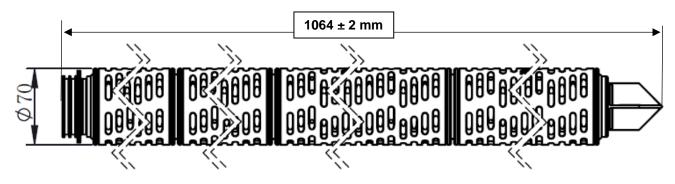

MCS – 40 inches

Part Number	Internal Filter Length	End Modification	Gasket Material		
(inches)					
MCS2074S 40"		Code 7: 226 O-Ring with Spear	Silicone		
Find the complete list of available standard inlet and outlet fittings in page 8.					

Typical Filtered Volume:

Filter Size	Filter Area	Volume to be Filtered (Thru-Put in Liters)	
5"	$0.35 m^2$	50 L-200 L	
10"	$0.7 m^2$	100 L-400 L	
20"	$1.4 m^2$	200 L-800 L	
30"	2.8 m ²	400 L-1600 L	
40"	5.6 m ²	800 L-3200 L	

*Volume Filtered is an estimated range for Sterile compounding, Pharmaceutical & Biological applications. The Filtration Volume is significantly influenced by your solution's viscosity, active ingredients, etc., properties. The best way to determine the exact filtration volume per filter is to try it with your solution.

Filter Integrity:

Revolutionary systems for your industry

The finished product was sampled and shown to exhibit a minimum bubble point of ≥18 psi (1.2 bar) in 60%IPA / 40% water and \geq 50 psi (3.5 bar) in DI Water (at 22°C).





International Filter Products

Revolutionary systems for your industry

Sterilization:

Autoclave: The filters can be sterilized by autoclaving for up to 50 cycles at 125 °C (257 °F) for 30 minutes.

Steam-In-Place: The filters can also be sterilized by steam-in-place procedure up to 30 cycles at 135 °C (275 °F) for 30 minutes at less than 0.3 bar differential pressure.

Sanitization: The filters can also be sanitized by hot water or common chemicals that are compatible with filter components.

Regulatory Compliance 21CFR Part 177 & USP <88>:

The filters are constructed with polypropylene resins and filtration media in compliance with 21CFR Part 177 of the US Code of Federal Regulations and USP Class VI Biological USP <88>Test for Plastic.

USP <85> Bacterial Endotoxins:

The filters were tested to confirm that an aqueous extraction of this product contains <0.25 EU/ml as determined by the Limulus Amebocyte Lysate (LAL) Test

USP <645> Conductivity USP:

Effluent is tested during the manufacturing process and shown to meet the requirements for USP Sterile Water for Injection for conductivity.

Bacterial Retention ASTM F838-05:

The filters are quantitatively retentive towards *Brevundimonas diminuta* (ATCC #19146) at a minimum challenge level of 10^7 CFU per cm^2 of filtration area, consistent with ASTM F838-05

USP <87> / ISO 10993-5 Cytotoxicity:

Extract from this product is non-Cytotoxic

Hemolysis ASTM F756-17:

Extract from this product is non-Hemolytic

Human and Veterinarian Use:

Revolutionary systems for your industry

The product is safe for Human and Veterinarian use. CGMP CFR part 210 & 211, additional requirement of 21 CFR part 600 and 21 CFR part 680 are applicable to the aseptic manufacturing process.

Animal-Derived Components & TSE/BSE Risk:

International Filter Products

No animal-derived material is intentionally added or used during the manufacture of this product.

Shelf Life:

The MCS capsules have a shelf life of 3 years from the date of manufacture.



FP International Filter Products Revolutionary systems for your industry

MCS - Series PES Cartridge

(Pharmaceutical Grade)

Ordering Guide

M C S [1] [2] [3] [4] 1 [5] [6] -PH [7]						
[1]	[2]	[3]	[4]	[5]	[6]	[7]
Micron Rating	End Modification	Length	O-Ring/Gasket Material	Optional	Pre-Filter	Sterilization
 04 : 0.04 µm 10 : 0.10 µm 20 : 0.20 µm 50 : 0.45 µm 65 : 0.65 µm 80 : 0.80 µm 1X : 1.20 µm Note: All final membranes are sterilizing grade.	 0: 222 O-Ring Flat 5: 222 O-Ring Spear 6: 226 O-Ring Flat 7: 226 O-Ring Spear F: DOE Flat Gasket S: SOE Flat Gasket 	 1: 10-Inch 2: 20-Inch 3: 30-Inch 4: 40-Inch 5: 5-Inch 	 S: Silicone E: EPDM N: Buna-N P: Peroxide Cured EPDM Q: Platinum Cured Silicon T: TEV or FEP Gasket V: Viton 	 -5 : SS Insert -RI : RFID Chip 	 DS065 : Highly Asymmetric 0.65µm DS045 : Highly Asymmetric 0.45µm S080 : 0.8µm 	 -ETO: Ethylene Oxide Sterilization Note: Only available for 5-in & 10-in Cartridges

Example:

MCS2071S-PH-ETO

PureFlo MCS Series Cartridge, 0.2µm PES membrane (Sterilizing Grade), 10.0-inch Long, Filtration Area 0.7 m2, 226 O-Rings with Spear, Silicone Double O-Rings, Stainless Steel Insert, ETO Sterilized (Factory Sterilized)

Part Number

Description





Special Configuration:

Layer Option:

- Single Layer
- Dual Layer

Micron Rating Option:

- 0.04 Micron
- 0.1 Micron
- 0.2 Micron
- 0.45 Micron
- 0.65 Micron
- 0.8 Micron
- 1.2 Micron

End Modification Option:

- 0:222 O-Ring Flat
- 3: 222 O-Ring w /tabs Spear
- 5:222 O-Ring Spear
- 6:226 O-Ring Flat
- 7:226 O-Ring Spear
- 8: 223 O-Ring Flat
- F : DOE Flat Gasket
- S : SOE Flat Gasket

O-Ring/Gasket Material Option:

- S : Silicone
- E : EPDM
- N : Buna-N
- P : Peroxide Cured EPDM
- Q : Platinum Cured Silicon
- T: TEV or FEP Gasket
- V:Viton

FP International Filter Products Revolutionary systems for your industry